O₂ ISM Qualification Kit

Instruction manual Bedienungsanleitung Instructions d'utilisation

O₂ ISM Qualification Kit
52 300 440 F

METTLER TOLEDO

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O₂ ISM Qualification Kit

Instruction manual

ISM and InPro are trademarks of the METTLER TOLEDO Group.

Contents

	Page
1 Safety	5
2 Important notes	6
2.1 Notes on operating instructions	6
2.2 Intended use	6
2.3 Safety instructions	7
2.4 Environmental protection	9
2.5 Examples of some typical applications	9
2.6 Use in Ex-zones	10
2.7 Ex-classification ATEX	11
2.7.1 Introduction	11
2.7.2 Rated data	11
2.7.3 Special conditions	12
2.8 Ex-classification FM approved	13
2.8.1 Introduction	13
2.8.2 Rated data	14
2.8.3 Special conditions	14
2.8.4 Applied Standards	15
3 Product Description	16
4 Operation	17
4.1 Connecting the transmitter to the O ₂ ISM Qualification Kit	17
5 Product Specification	20
5.1 Simulator, Pre-registered Values	20
5.2 Sensor Diagnostics Data	21
5.3 Warnings and Alarms	23
6 Installation	25
6.1 Dimensions	25
7 Ordering Information	26
8 Protocol	27

1 Safety

This instruction manual lists the main points of how to use the **O₂ ISM™ Qualification Kit**, from METTLER TOLEDO for best results and in the intended manner.

This **instruction manual**, and the **safety requirements** in particular, **must be observed** by the **operator and all other persons** working with this equipment.

The instruction manual must always be stored close at hand, in a place accessible to all people working with an ISM O₂ sensor and the O₂ ISM Qualification Kit.

Warning notices and symbols

This instruction manual identifies safety instructions and additional information by means of the following symbols:

This symbol draws attention to safety instructions and **warnings of potential danger** which, if neglected, could result in injury to persons and/or damage to property.

This symbol identifies **additional information and instructions** which, if neglected, could lead to defects, inefficient operation and possible loss of product.

2 Important notes

2.1 Notes on operating instructions

These operating instructions contain all the information needed for safe and proper use of the O₂ Qualification Kit.

The operating instructions are intended for personnel entrusted with the operation and maintenance of oxygen sensors. It is assumed that these persons are familiar with the equipment in which the sensors are installed.

Warning notices and symbols

This instruction manual identifies safety instructions and additional information by means of the following symbols:

This symbol draws attention to **safety instructions and warnings of potential danger** which, if neglected, could result in injury to persons and/or damage to property.

This symbol identifies **additional information and instructions** which, if neglected, could lead to defects, inefficient operation and possible loss of production.

2.2 Intended use

The METTLER TOLEDO O₂ ISM Qualification Kit is intended solely for the qualification of O₂ measuring systems as described in this instruction manual.

Any use of this kit which differs from or exceeds the scope of use described in this instruction manual will be regarded as inappropriate and incompatible with the intended purpose.

The manufacturer/supplier accepts no responsibility whatsoever for any damage resulting from such improper use. The risk is borne entirely by the user/operator.

Other prerequisites for appropriate use include:

- Compliance with the instructions, notes and requirements set out in this instruction manual.
- Acceptance of responsibility for regular inspection, maintenance and functional testing of all associated components, also including compliance with local operational and plant safety regulations.
- Compliance with all information and warnings given in the documentation relating to the products used in conjunction with the sensor (housings, transmitters, etc.).
- Observance of all safety regulations governing the equipment in which the sensor is installed.
- Correct equipment operation in conformance with the prescribed environmental and operational conditions, and admissible installation positions.
- Consultation with METTLER TOLEDO Process Analytics in the event of any uncertainties.

2.3 Safety instructions

- The plant operator must be fully aware of the potential risks and hazards attached to operation of the particular process or plant. The operator is responsible for correct training of the workforce, for signs and markings indicating sources of possible danger, and for the selection of appropriate, state-of-the-art instrumentation.
- It is essential that personnel involved in the commissioning, operation or maintenance of these simulators or of any of the associated equipment (e.g. housings, transmitters, etc.) be properly trained in the process itself, as well as in the use and handling of the associated equipment. This includes having read and understood this instruction manual.

- The safety of personnel as well as of the plant itself is ultimately the responsibility of the plant operator. This applies in particular in the case of plants operating in hazardous zones.
- The oxygen sensor simulators and associated components have no effect on the process itself and cannot influence it in the sense of any form of control system.
- Where specific safeguards such as locks, labels, or redundant measuring systems are necessary, these must be provided by the plant operator.
- A defective simulator must neither be installed nor put into service.
- No modifications to the simulators and the accessories are allowed. The manufacturer accepts no responsibility for damage caused by unauthorized modifications. The risk is borne entirely by the user.
- When using cables for simulator connection not supplied and recommended by the manufacturer the following standards need to be considered:
 - US National Electrical Code® (ANSI/NFPA 70 [NEC®]), where applicable.
 - Canadian Electrical (CE) Code® (CEC Part 1, CAN/CSA-C22.1), where applicable.
- For guidance on US installations, see ANSI/ISA-RP12.06.01, Installation of Intrinsically Safe Systems for Hazardous (Classified) Locations.
- Control room equipment connected to intrinsically safe associated apparatus in the US should not use or generate more than the specified Um of the associated apparatus.
- Care must be taken during installation to avoid impacts or friction that could create an ignition source.
- Tampering and replacement with non-factory components may adversely affect the safe use of the system.

- Insertion or withdrawal of removable electrical connectors or modules is to be accomplished only when the area is known to be free of flammable vapors.
- **WARNING!**
INTRINSICALLY SAFE APPARATUS CAN BE A SOURCE OF IGNITION IF INTERNAL SPACINGS ARE SHORTED OR CONNECTIONS OPENED.
- **WARNING!**
DO NOT OPEN WHEN AN EXPLOSIVE ATMOSPHERE IS PRESENT.
- **WARNING!**
SUBSTITUTION OF COMPONENTS MAY IMPAIR INTRINSIC SAFETY.
- **WARNING!**
SUBSTITUTION OF COMPONENTS MAY IMPAIR SUITABILITY OF THE EQUIPMENT.
- **WARNING!**
FOR CONNECTION ONLY TO NON-FLAMMABLE PROCESSES.
- **WARNING!**
TO MAINTAIN THE ENCLOSURE IP66 RATING, THE CONNECTOR MUST BE FULLY ENGAGED.

2.4 Environmental protection

Waste electrical products should not be disposed of with household waste. Please recycle where facilities exist. Check with your Local Authority or retailer for recycling advice.

2.5 Examples of some typical applications

Below is a list of examples of typical fields of application for the oxygen sensor simulators. This list is not exhaustive:

- Tool for qualification of O₂ measuring systems
- Service tool for quick checks and loop troubleshooting

2.6 Use in Ex-zones

Attention!

For an installation in Ex-zones please read the guidelines following hereafter:

Ex-classification ATEX

⊕ II 1/2G Ex ia IIC T6...T3 Ga/Gb

⊕ II 1/2D Ex ia IIIC T26 °C...T161 °C Da/Db

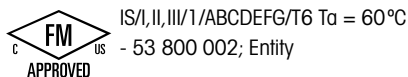
Number of the test certificate:

SEV 14 ATEX 0169X

IECEX SEV 14.0026X

CML 22 UKEX 2212X

Ex-classification FM approved:

Project ID: 3021227

FM certificate number: FM16US0034X
FM18CA0021X

2.7 Ex-classification ATEX

2.7.1 Introduction

According to Directive 2014/34/EU¹⁾ (ATEX 114) Appendix I, oxygen ISM sensor simulators are devices group II, category 1/2G and according to RL 99/92/EG²⁾ (ATEX 137) may be used in zones 0/1 or 1/2 or 1 or 2 and gas groups IIA, IIB and IIC that are potentially explosive due to combustible substances in above stated temperature classes.

For use / installation, the requirements of EN 60079-14 must be observed.

According to Directive 2014/34/EU¹⁾ (ATEX 114) Appendix I, oxygen ISM sensor simulators are devices group II, category 1/2D and according to RL 99/92/EG²⁾ (ATEX 137) may also be used in zones 20/21 or 21 or 22 that contain combustible dusts.

The digital sensor circuit is part of a common intrinsically safe system and is for operation connected to a separately certified transmitter.

The digital sensor circuit as part of an intrinsically safe system is isolated from the not-intrinsically safe electric circuits up to a maximum rated voltage of 375 V and from grounded parts up to a maximum rated voltage of 30 V.

2.7.2 Rated data

Measuring circuit:

Intrinsically safe ignition protection class Ex ia IIC only for connection to a certified intrinsically safe circuit.

Digital O₂ oxygen sensor simulator maximum values:

Two-wire current circuit: $U_i \leq 16 \text{ V}$ $L_i = \text{negligible}$
 $I_i \leq 30 \text{ mA}$ $C_i = \text{negligible}$
 $P_i \leq 50 \text{ mW}$

¹⁾ For UK statutory Requirements SI 2016 No. 1107

²⁾ For UK regulations 7 and 11 of the Dangerous Substances and Explosive Atmospheres Regulations 2002 (DSEAR).

Note: The above maximum values are each the total of all individual circuits of the associated intrinsically safe power supply and transmitter.

2.7.3 Special conditions

- The relationship between the maximum permissible ambient or media temperature and temperature class, for category 1G applications, zone 0, is shown in the following table:

Temperature class	Max. environment resp. media temperature
T 6	68 °C
T 5	80 °C
T 4	108 °C
T 3	160 °C

- The relationship between the maximum permissible ambient or media temperature and temperature class, for category 1G applications, zone 0, is shown in the following table:

Temperature class	Max. environment resp. media temperature
T69°C	68 °C
T81°C	80 °C
T109°C	108 °C
T161°C	160 °C

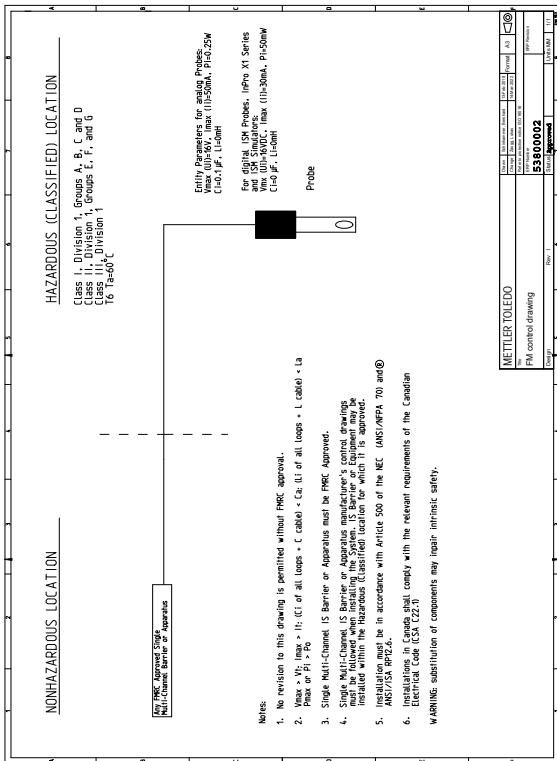
- The capacitance and inductance of the connecting cable has to be considered.

2.8 Ex-classification FM approved

2.8.1 Introduction

The following FM control drawing and the standards listed in section 2.7.4 must be observed, where applicable:

2.8.2 Rated data

For rated data please observe section 2.7.2.

2.8.3 Special conditions

In type of protection intrinsically safe apparatus, the analytical probes (oxygen ISM sensor simulators) equipment is designated with the following specific conditions of use.

1. The oxygen ISM sensor simulators shall be installed in compliance with the enclosure, mounting, spacing and segregation requirements of the ultimate application, including a tool removable cover.
2. Process temperature no greater than +130 °C.
3. Potential Electrostatic Charging Hazard – To prevent the risk of electrostatic sparking, the non-metallic surface should only be cleaned with a damp cloth.

2.8.4 Applied Standards

United States Standards

Titel	Number	Issue Date
Approval Standard for Electrical Equipment for Use in Hazardous (Classified) Locations – General Requirements.	FM Class 3600	2022
Approval Standard for Intrinsically Safe Apparatus and Associated Apparatus for Use in Class I, II & III, Division 1, Hazardous (Classified) Locations.	FM Class 3610	2021
Approval Standard for Electrical Equipment for Measurement, Control and Laboratory Use.	FM Class 3810	2021
Explosive Atmospheres – Part 0: Equipment – General Requirements.	ANSI/UL 60079-0	2022
Explosive Atmospheres – Part 11: Equipment Protection by Intrinsic Safety “i”.	ANSI/UL 60079-11	2018
Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use – Part 1: General Requirements.	ANSI/UL 61010-1	2018

Canadian Standards

Titel	Number	Issue Date
Explosive Atmospheres – Part 0: Equipment – General Requirements.	CAN/CSA-C22.2 No. 60079-0	2019
Explosive Atmospheres – Part 11: Equipment Protection by Intrinsic Safety “i”.	CAN/CSA-C22.2 No. 60079-11	R2018
Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use – Part 1: General Requirements.	CAN/CSA-C22.2 No. 61010-1	R2017

3 Product Description

The O₂ ISM Qualification Kit is a set of 5 different sensor plug-in heads and allows the simulation of the following sensor types with predefined values and errors (not changeable by the user) to assure safe use as qualification tool:

- InPro™ 6850 i
- InPro 6900i
- InPro 6950 i

Each sensor plug-in head corresponds to one of the mentioned METTLER TOLEDO sensors and delivers the complete data of an ISM sensor.

The O₂ ISM Qualification Kit has basically 2 functions:

- Tool for the qualification of measuring systems.
- Service tool for quick checks and loop trouble-shooting.

The data indicated (primary/ secondary) and transmitted are as follow:

- Main measured value
- Temperature
- Error message according to indication.

All standard process and wear data of the sensor are also available in each plug-in head and are also sent to the transmitter or the PLC as well as to the iSense software.

4 Operation

4.1 Connecting the transmitter to the O₂ ISM Qualification Kit

To qualify your measuring systems proceed as follows:

Please refer to the transmitter instruction manual for the exact description of the transmitter operation.

Step 1: Disconnect the ISM sensor from the measuring point.

Step 2: Connect one of the O₂ ISM Qualification Kit plug-in heads to the transmitter via the standard connecting cable.

Step 3: The preregistered values then are displayed on the transmitter.

Step 4: Check if the displayed values by the transmitter or the PLC correspond to the predefined fixed values of the O₂ ISM Qualification Kit plug-in head.

See certification for exact values and tolerances.

Step 5: In case you are using the Error simulators, so check that the appropriate errors are correctly displayed by the transmitter or the PLC.

Step 6: Report results into the attached spreadsheet "Protocol for O₂ ISM Qualification Kit".

Please note: the O₂ ISM Qualification Kit simulates the behavior of a sensor at several temperatures. All commonly used tables programmed in each transmitter for solubility, salinity correction, etc. apply standardized values, e.g. temperature 25 °C (77 °F), pressure 1.013 bar (14.69 psi) and a relative humidity 100 %. The transmitter calculates back the value that the sensor would deliver under these reference conditions. This compensated value may differ from the direct current delivered by the simulator if for example the calibration pressure (manually tipped or automatically measured by some transmitter) differs from 1.013 bar (14.69 psi) or if the relative humidity typed in is different from 100 %.

Problem solving: If any of the tests described above failed, you should change the connecting cable and repeat the test. If the test still shows the same result, the transmitter or the cable is defective. If you are using a transmitter with current outputs, so ensure the outputs are correctly wired.

You must **never do a calibration** (the simulators are read-only).

You must **never increment autoclave cycles** (always set "NO" to cycle counter), this is only needed if autoclave has been set to a value higher than 0.

Different transmitters may display different values and errors.

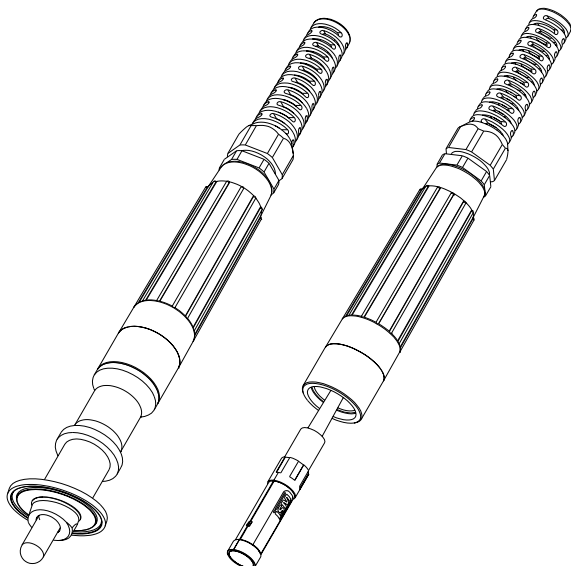
Out of range behavior is also different depending on the transmitter in use thus **different values/readings may occur.**

Set the transmitter input filter to off (only for SIM O₂ TGL, SIM O₂ ppb TGL and SIM O₂ Trace TGL).

Do not use the O₂ ISM plug-in heads after the expiry date shown on the label of each plug-in head. Please send them back to METTLER TOLEDO for recertification.

Measuring point with
METTLER TOLEDO Electrode

Connecting O₂ ISM
Qualifying plug-in head

Environment operating and storing conditions:

Permissible ambient temperature (operation):

- 25 °C to 60 °C (- 13 °F to 140 °F)

Permissible ambient temperature (storage/transport):

- 25 °C to 60 °C (- 13 °F to 140 °F)

5 Product Specification

5.1 Simulator, Pre-registered Values

Order No.	Product Designation	Legend	OXY Value	Temp.
Standard (InPro 6850i) SIM O ₂ Kit			Order No. 52 300 416	
52 300 417	SIM O ₂ ZERO	SIM O ₂ ZERO	0%	25 °C
52 300 418	SIM O ₂ AIR	SIM O ₂ AIR	20.95 Vol% 100% (air) sat	40 °C
52 300 419	SIM O ₂ TGL	Toggle Mode 30 sec	0% & 100%	25 °C & 40 °C
52 300 420	SIM O ₂ ERR1	9999 % ...% sat	9999 % ...% sat	25 °C
52 300 421	SIM O ₂ ERR2	9999 % ...% sat	83 % ...% sat	90 °C
Standard (InPro 6900i) SIM O ₂ ppb Kit			Order No. 52 300 422	
52 300 423	SIM O ₂ ppb ZERO	SIM O ₂ ppb ZERO	0%	5 °C
52 300 424	SIM O ₂ ppb AIR	SIM O ₂ ppb AIR	20.95 Vol% 100% (air) sat	25 °C
52 300 425	SIM O ₂ ppb TGL	Toggle Mode 30 sec	0% & 100%	5 °C & 25 °C
52 300 426	SIM O ₂ ppb ERR1	9999 % ...% sat	9999 % ...% sat	25 °C
52 300 427	SIM O ₂ ppb ERR2	9999 % ...% sat	83 % ...% sat	90 °C
Standard (InPro 6950i) SIM O ₂ Trace Kit			Order No. 52 300 428	
52 300 429	SIM O ₂ Trace ZERO	SIM O ₂ Trace ZERO	0%	5 °C
52 300 430	SIM O ₂ Trace AIR	SIM O ₂ Trace AIR	20.95 Vol% 100% (air) sat	25 °C
52 300 431	SIM O ₂ Trace TGL	Toggle Mode 30 sec	0% & 100%	5 °C & 25 °C
52 300 432	SIM O ₂ Trace ERR1	9999 % ...% sat	9999 % ...% sat	25 °C
52 300 433	SIM O ₂ Trace ERR2	9999 % ...% sat	83 % ...% sat	90 °C

5.2 Sensor Diagnostics Data

All O₂ SIM's except 52 300 421 (SIM O₂ ERR2), 52 300 427 (SIM O₂ PPB ERR2), 52 300 433 (SIM O₂ TRACE ERR2)

Sensor Diagnostics O₂	Transmitters M400 / M420 / M800
Time to maintenance (TTM)	100 d
Life time indicator (DLI)	2400 d
Adaptive Cal Timer (ACT)	50 d
Sensor operating time	300 d
Autoclaving cycles	1
CIP cycles	3
SIP cycles	2
Max temperature	130 / 121 °C dd.mm.yyyy (date of production)

For SIM 52 300 421 (SIM O₂ ERR2), 52 300 427 (SIM O₂ PPB ERR2), 52 300 433 (SIM O₂ TRACE ERR2)

Sensor Diagnostics O₂	Transmitters M400 / M420 / M800
Time to maintenance (TTM)	0 d
Life time indicator (DLI)	2300 d
Adaptive Cal Timer (ACT)	0 d
Sensor operating time	300 d
Autoclaving cycles	1
CIP cycles	3
SIP cycles	2
Max temperature	130 / 121 °C dd.mm.yyyy (date of production)

All O₂ SIM except 52 300 421 (SIM O₂ ERR2), 52 300 427 (SIM O₂ PPB ERR2), 52 300 433 (SIM O₂ TRACE ERR2)

Sensor Wear Monitor O₂	Transmitter M700 % bar
Membrane body wear	~100 %
Inner body wear	~10 %
Operating time	300 d
Network diagram sensor wear	50 %

For SIM 52 300 421 (SIM O₂ ERR2), 52 300 427 (SIM O₂ PPB ERR2), 52 300 433 (SIM O₂ TRACE ERR2)

Sensor Wear Monitor O₂	Transmitter M700 % bar
Membrane body wear	~100 %
Inner body wear	~15 %
Operating time	300 d
Network diagram sensor wear	25 %

5.3 Warnings and Alarms

Warnings and Alarms	Transmitters M300 / M400 / M420
52 300 420 (SIM O ₂ ERR1)	Error O ₂ slope > – 30 nA
52 300 426 (SIM O ₂ ppb ERR1)	Error O ₂ slope > – 220 nA
52 300 432 (SIM O ₂ Trace ERR1)	Error O ₂ slope > – 2000 nA
52 300 421 (SIM O ₂ ERR2)	Calibration required Maintenance required
52 300 427 (SIM O ₂ ppb ERR2)	Calibration required Maintenance required
52 300 433 (SIM O ₂ Trace ERR2)	Calibration required Maintenance required

Warnings and Alarms	Transmitters M700
52 300 420 (SIM O ₂ ERR1)	Fail Measurement Range Warning Slope
52 300 426 (SIM O ₂ ppb ERR1)	Fail Measurement Range Warning Slope
52 300 432 (SIM O ₂ Trace ERR1)	Fail Measurement Range Warning Slope
52 300 421 (SIM O ₂ ERR2)	Fail Measurement Range Warning Sensor wear Warning Adjustment required
52 300 427 (SIM O ₂ ppb ERR2)	Fail Measurement Range Warning Sensor wear Warning Adjustment required
52 300 433 (SIM O ₂ Trace ERR2)	Fail Measurement Range Warning Sensor wear Warning Adjustment required


Warnings and Alarms	Transmitters M800
52 300 420 (SIM O ₂ ERR1)	Error O ₂ hi Slope > – 30 nA
52 300 426 (SIM O ₂ ppb ERR1)	Error O ₂ lo Slope > – 220 nA
52 300 432 (SIM O ₂ Trace ERR1)	Error O ₂ Trace Slope > – 2000 nA
52 300 421 (SIM O ₂ ERR2)	Maint Required Cal Required
52 300 427 (SIM O ₂ ppb ERR2)	Maint Required Cal Required
52 300 433 (SIM O ₂ Trace ERR2)	Maint Required Cal Required

Please refer to your transmitter instruction manual for specific instructions.

6 Installation

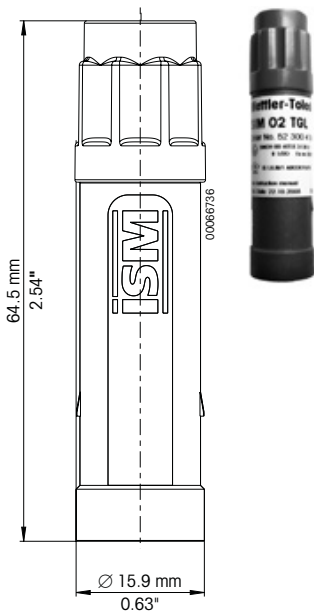6.1 Dimensions

Markings

The markings labeled on each pH ISM plugin head state the following information:


Simulator O2 ZERO		METTLER TOLEDO www.mt.com/gro	
P/N:	52300417		Mettler-Toledo GmbH Im Hackacker 15 CH-8902 Urdorf Switzerland
S/N:	1234567		SEV 14 ATEX 0169X IECEx SEV 14.0026X CML 22 UKEX 2212X See instruction manual!
	HAZ LOC, IS/II/III/1/ ABCDEF/G/T6; Ta<+60°C FM16US0034X FM18CA0021X		
Cal. Date:	02.09.2022		1258 UK CA 2503
Cal. Due:	01.09.2023		
HW: 02.00.00	FW: SIM-05.00.00		

7 Ordering Information

Order No.	Designation	Order No.	Designation
52 300 410	SIM pH Kit	52 300 428	SIM O₂ Trace Kit
52 300 411	SIM pH4	52 300 429	SIM O ₂ Trace ZERO
52 300 412	SIM pH7	52 300 430	SIM O ₂ Trace AIR
52 300 413	SIM pH TGL	52 300 431	SIM O ₂ Trace TGL
52 300 414	SIM pH ERR1	52 300 432	SIM O ₂ Trace ERR1
52 300 415	SIM pH ERR2	52 300 433	SIM O ₂ Trace ERR2
52 300 416	SIM O₂ Kit	52 300 422	SIM O₂ ppb Kit
52 300 417	SIM O ₂ ZERO	52 300 423	SIM O ₂ ppb ZERO
52 300 418	SIM O ₂ AIR	52 300 424	SIM O ₂ ppb AIR
52 300 419	SIM O ₂ TGL	52 300 425	SIM O ₂ ppb TGL
52 300 420	SIM O ₂ ERR1	52 300 426	SIM O ₂ ppb ERR1
52 300 421	SIM O ₂ ERR2	52 300 427	SIM O ₂ ppb ERR2
30 031 035	SIM CO₂ Kit		
30 034 395	SIM CO ₂ 15 mbar		
30 038 175	SIM CO ₂ 950 mbar		
30 038 973	SIM CO ₂ TGL		
30 038 974	SIM CO ₂ ERR1		
30 038 975	SIM CO ₂ ERR2		

The single simulators are not available for sale.

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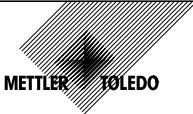A copy of this form in A4 format is delivered with the O₂ ISM Qualification Kit.

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2.7.4 Angewandte Normen

US-Normen

Titel	Nummer	Ausgabe Datum
Approval Standard for Electrical Equipment for Use in Hazardous (Classified) Locations – General Requirements.	FM Class 3600	2022
Approval Standard for Intrinsically Safe Apparatus and Associated Apparatus for Use in Class I, II & III, Division 1, Hazardous (Classified) Locations.	FM Class 3610	2021
Approval Standard for Electrical Equipment for Measurement, Control and Laboratory Use.	FM Class 3810	2021
Explosive Atmospheres – Part 0: Equipment – General Requirements.	ANSI/UL 60079-0	2022
Explosive Atmospheres – Part 11: Equipment Protection by Intrinsic Safety “I”.	ANSI/UL 60079-11	2018
Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use – Part 1: General Requirements.	ANSI/UL 61010-1	2018

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Titel	Nummer	Ausgabe Datum
Explosive Atmospheres – Part 0: Equipment – General Requirements.	CAN/CSA-C22.2 No. 60079-0	2019
Explosive Atmospheres – Part 11: Equipment Protection by Intrinsic Safety “I”.	CAN/CSA-C22.2 No. 60079-11	R2018
Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use – Part 1: General Requirements.	CAN/CSA-C22.2 No. 61010-1	R2017

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Titre	Numéro	Date de publication
Approval Standard for Electrical Equipment for Use in Hazardous (Classified) Locations – General Requirements.	FM Class 3600	2022
Approval Standard for Intrinsically Safe Apparatus and Associated Apparatus for Use in Class I, II & III, Division 1, Hazardous (Classified) Locations.	FM Class 3610	2021
Approval Standard for Electrical Equipment for Measurement, Control and Laboratory Use.	FM Class 3810	2021
Explosive Atmospheres – Part 0: Equipment – General Requirements.	ANSI/UL 60079-0	2022
Explosive Atmospheres – Part 11: Equipment Protection by Intrinsic Safety "i".	ANSI/UL 60079-11	2018
Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use – Part 1: General Requirements.	ANSI/UL 61010-1	2018

Normes canadiennes

Titre	Numéro	Date de publication
Explosive Atmospheres – Part 0: Equipment – General Requirements.	CAN/CSA-C22.2 No. 60079-0	2019
Explosive Atmospheres – Part 11: Equipment Protection by Intrinsic Safety "i".	CAN/CSA-C22.2 No. 60079-11	R2018
Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use – Part 1: General Requirements.	CAN/CSA-C22.2 No. 61010-1	R2017

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All rights reserved
Protocole pour set de qualification ISM O₂ © Mettler-Toledo AG, MarCom, Im Hackacker 15, CH-8902 Urdorf. 03/2009 1